S*PURE Application Note - Environmental

Triazine Herbicides in Water On SEClute SPE Column



SPE Conditions

Procedure using SEClute™ C18-Fast columns, 500mg/3mL, (Part #5138758)

Sample Treatment: Spiked 500mL of water spiked with 10ug each triazine herbicide (simazine, atrazine, propazine, terbuthylazine, and cyanazine).

Conditioning: 2 x 2mL acetone, 2 x 2mL methanol, 2 x 2mL water. Column must not run dry!

Sample Application: Attached a 75mL reservoir (Part #210575) to the SPE tube with a syringe adapter (Part # 210705) and aspirate the sample through the tube.

Elution: Elute triazine herbicides with 3 x 1mL acetone. Adjust sample volume to 1mL by evaporating off acetone with a warm nitrogen stream.

GC Conditions

Injected 1.6uL of above sample onto AT™ 5ms column, 30m x 0.25mm i.d., 0.25µm film thickness (Part #15807) under the following conditions:

Oven: 80°C (0 minutes)-280°C (0 minutes) at 10°C/min


Injector: 250°C

FID: 310°C, 14.5psig helium

Linear Velocity: 29cm/sec Flow: 1.1mL/minute

Split Ratio: 50:1

SVF: 55mL/minute (all per Agilent 6890)

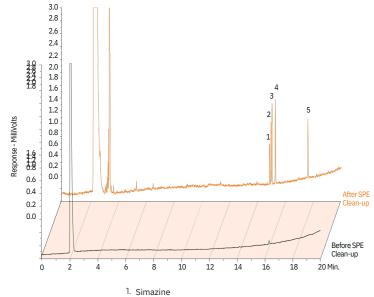

2. Atrazine 3. Propazine 4. Terbuthylazine 5. Cyanazine

www.sstarpure.com

AT" is a trademark of Altech Associates, Inc. GRACE* and GRACE DAVISON* are trademarks, registered in the United States and/or other countries, of W. R. Grace & Co.-Conn, SEClute** is a trademark of S*Pure Pte Ltd, Singapore. This trademark list has been compiled using available published rformation as of the publication date of this brochure and may not accurately reflect current trademark. An enership, All rights reserved. The information presented herein is derived from our testing and experience. It is direred for your consideration and verification. Since operating conditions vary isofficative and are on under our control wardiscipant in the publication is a trademark of the control of the contr

S*Pure Pte Ltd

55 Toh Guan Road East #03-05 Unitech Centre Singapore 608601 Tel: +65 6563 3901 Email: corporate@sstarpure.com 2024-08, M212