S*PURE Application Note - Environmental

Chorlinated Pesticides in Vegetables On SEClute SPE Column



SPE Conditions

Procedure using SEClute[™] C18-Max SPE tube, 500mg/6mL (Part #5138767)

Sample Treatment: Homogenize 5 grams green bell pepper with 25mL methanol, spike with 0.5 μ g each of 16 chlorinated pesticides. Filter with 75mL filter column (Part #210775). Dilute to 200mL with DI water and shake.

Conditioning: 2 x 5mL methanol, then 5mL DI water.

Sample Application: Attach 75mL reservoir (Part #210575) with a syringe adapter (Part #210705) and apply entire sample.

Elution: Elute chlorinated pesticides with 2 x 1mL n-hexane/CH2Cl2 (70:30), evaporated this to dryness and reconstituted with 1mL n-hexane.

GC Conditions

Capillary column (Part #15807), AT[™] 5ms, 30m x 0.25mm i.d., 0.25µm film thickness.

Oven: 80°C (1min) to 100°C at 15°C/min, to 160°C at 10°C/min, to 275°C at 5°C/min, column head pressure at 14.0psi on the EPC-1000, dead time of 1.91 minutes Linear Velocity: 26cm/sec


Flow: 0.77mL/minute

Split Vent Flow: 50mL/minute

Split Ratio: 65:1

Injector: 250°C

ECD: 310

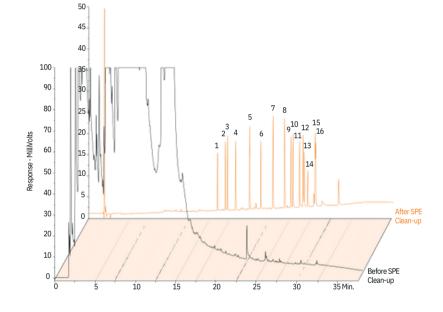

1. a-BHC
2.b-BHC
3.g-BHC
4.d-BHC
5.Heptachlor
6.Aldrin
7.Heptachlor Epoxide
8.Endosulfan I

9. p,p'-DDE 10. Dieldrin 11.Endrin 12.Endosulfan II 13.p,p'-DDD 14.Endrin Aldehyde 15.p,p'-DDT 16.Endosulfan Sulfate

www.sstarpure.com

AT^{**} is a trademark of Alltech Associates, Inc. GRACE[®] and GRACE DAVISON[®] are trademarks, registered in the United States and/or other countries, of W. R. Grace & Co.-Conn. SEClute[™] is a trademark of S^{*}Pure Pte Ltd, Singapore. This trademark is thas been compiled using available published information as of the publication date of this brochure and may not accurately reflect current trademark ownership. All highs reserved. The information presented herein is derived from our testing and experience. It is offered for your consideration and verification. Since operating conditions vary significantly, and are not under our control, we disclarate may be obtained from the use of our products. S^{*}Pure reverse the right to charge prices and/or specifications. Which up ion onflication.

S*Pure Pte Ltd

55 Toh Guan Road East #03-05 Unitech Centre Singapore 608601 Tel: +65 6563 3901 Email: corporate@sstarpure.com 2024-08, M206